

DHM's Haupt's Adhesive Concentrate

DHM's Haupt's Adhesive Concentrate saves time, money, and the environment by eliminating prep work, clean up, and waste. Conveniently packaged in 30mL disposable polypropylene vials, your 15mL concentrated aliquot (gently heated to liquid form using a lab safe microwave, water bath, or oven) quickly turns into 30mL of working solution by simply adding 15mL of diluted EtOH (1:1 ratio w/ 50% EtOH recommended). After inverting several times (approximately 30 seconds recommended) to thoroughly mix the solution, your Haupt's Adhesive solution is now ready for use. It's that simple. Each 30mL vial of working solution "coats" 200-300 slides (depending on slide size and number of drops used per slide) and dries clear within 5-10 minutes. Additionally, use of this adhesive, in conjunction with DHM's Resin Section Slide Press, optimizes both section quality and section retention throughout even the harshest of staining procedures. DHM's Haupt's Adhesive Concentrate has a shelf-life of greater than 1 year (stored @ 2-8C) and the diluted working solution has a shelf-life of approximately 1 week (stored @ room temperature).